

Magnetic Separation System, 60-position



Description

A magnetic separation system designed to hold and separate to the side of the tubes magnetic particles in 12 X 75 mm tubes. The two-piece system consists of a test tube rack, which fits over a magnetic separation rack containing rare earth magnets. The two-piece design allows test tubes to be set up, incubated and magnetically separated without removing the assay tubes from the test tube holder. The unit accommodates up to 60 test tubes.

Use

Place test tubes in the test tube holder, place the test tube holder on top of the magnetic separation rack and allow separation to proceed from 2-10 minutes depending on the size of the particles. With the rack in place invert the unit to decant the tubes.

Application

Separation unit can be used for a variety of applications including cell sorting, purification of proteins, molecular biology, immunoassays, etc.

Warning

Strong magnetic field. The separator should be used and stored at a safe distance from magnetic media, videotapes, credit cards, etc.

Ordering Information

PN 472231 **Magnetic Separation System, 60-position**

Abraxis LLC

54 Steamwhistle Drive
Warminster, PA 18974
Phone: (215) 357-3911
FAX: (215) 357-5232
Email: info@abraxiskits.com
<http://www.abraxiskits.com>

